

Manutention :

Toute manutention de poids et masses étalons doit se faire en respectant des règles d'hygiène adaptées.

Éviter :











- de saisir les poids de classes E et F à mains nues
- d'exposer les poids à la poussière et aux salissures
- de stocker les poids en milieu humide
- de soumettre les poids à un fort champ magnétique
- de tirer ou de pousser les poids sur une surface



Retrouvez tous nos accessoires dans notre tarif



ACCESSOIRES POUR POIDS ET MASSES ÉTALONS DE CLASSES E & F

	<=50 g	100 à 200 g	500 g à 2 kg	5 à 50 kg	>50 kg
Pincette 10 cm - amagnétique					
Pincette 15 cm - amagnétique					
Fourches et poignées de préhension					
Gants					
Chiffon doux non pelucheux					
Kit de nettoyage de poids					



Entretien :

Aucun entretien est nécessaire.

ATTENTION : Toute intervention pouvant modifier la valeur de la masse est à proscrire, par exemple :

- polissage
- remise en peinture
- ouverture de la cavité d'ajustage

Stockage :

Il est conseillé de stocker les poids et masses étalons dans leurs coffrets d'origine. Les poids et masses étalons sans coffret sont à stocker dans des endroits appropriés.



Environnement	Stockage des poids de classes E & F	Stockage des poids de classe M
Température de stockage	20°C ± 5°C	20°C ± 10°C
Hygrométrie	50% ±25%	
Milieu	Hors poussière - Hors intempéries	

Handling :

Any handling of weights and masses shall be in accordance with adapted hygiene rules.

Avoid :








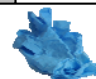


- taking weights of classes E and F with bare hands
- exposing the weights to dust and dirt
- storing weights in a humid atmosphere
- submitting weights to a strong magnetic fields
- pulling or pushing weights on a surface



Find all our accessories in our price-list



ACCESSORIES FOR STANDARD WEIGHTS AND MASSES OF CLASSES E AND F

	<=50 g	100 to 200 g	500 g to 2 kg	5 to 50 kg	>50 kg
Non-magnetic tweezer 10 cm					
Non-magnetic tweezer 15 cm					
Weight forks and handles					
Gloves					
Soft cloth					
Cleaning set					



Maintenance :

No maintenance is required.

Be careful : Any intervention that could change the value of the mass should be avoided, for example :

- polishing
- repainting
- opening the adjusting cavity

Storage :

It is recommended to store standard weights and masses in their original boxes. Weights and masses without box shall be stored in appropriate places.



Environment	Storage of weights of classes E and F	Storage of weights of class M
Storage temperature	20°C ± 5°C	20°C ± 10°C
Humidity	50% ±25%	
Dust	Off - Except bad weather	

F.AD.1401 2

ZWIEBEL SAS - SAINT JEAN SAVERNE - BP 50002 - F-67701 SAVERNE CEDEX

Tel. : +33 (0)3.88.71.53.10 - Fax: +33 (0)3.88.71.20.92

E-mail : commercial@zwibel.fr - www.zwibel.fr